

# Keysight Technologies

Topography and Recognition Imaging with PicoTREC,  
a Scanning Probe Microscopy Accessory for Mapping Target Molecules on  
a Sample Surface

Application Note

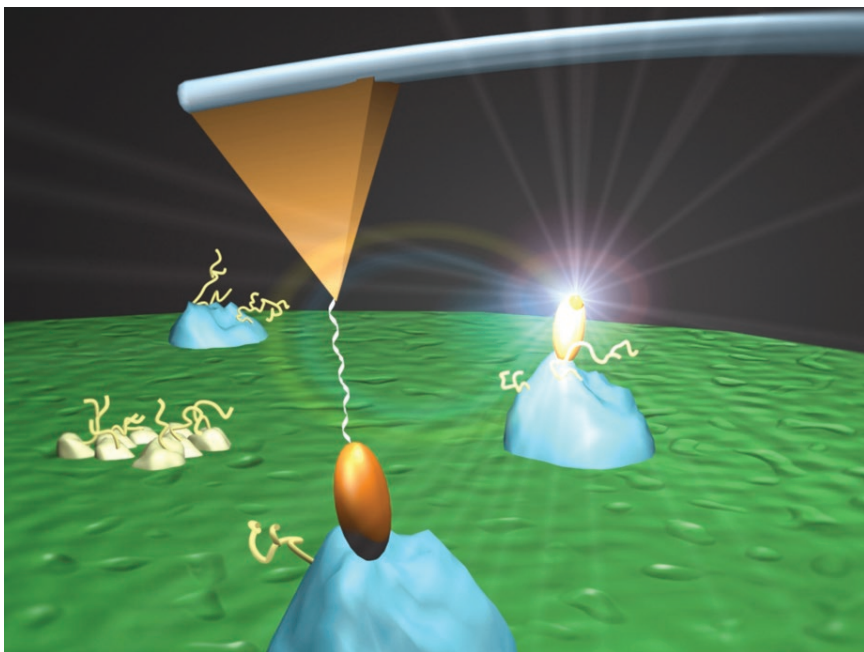


Figure 1. Schematic representation of an AFM tip measuring the nanomechanical properties of a flexible tether bound to a biological molecule (brown) recognizing its target (blue) on a surface. In TREC (topography and recognition) scanning mode, target molecules can be detected on the surface.

The development of the atomic force microscope (AFM) has allowed scientists and engineers to observe the details of molecular structures with unprecedented resolution and without the need for rigorous sample preparation or labeling. In addition to topographical imaging, AFM can sense nanomechanical and other fundamental properties of sample surfaces, including local adhesion and elasticity. Microscopic adhesion affects a variety of events from the behavior of ceramics and semiconductors to DNA replication and drug binding in the human body. The AFM offers a tool to study these important parameters on the nanometer scale using a technique that detects forces on the AFM probe tip as it approaches and retracts from a surface.

The Institute for Biophysics at the University of Linz, Austria, has developed a technique in which biological molecules are attached to the AFM tip on the end of a linker (Figure 1). The linker is an elastic PEG (polyethylene glycol) tether that gives the sensor molecule the freedom to reorient itself properly to bind to its target on the surface. When a recognition event occurs, the AFM detects the force required to break the bonds that are involved in molecular binding. This technique is being used to detect binding forces between antibody-antigen pairs, drug-receptors, and sense-antisense DNA complexes (1).

When combined with dynamic force microscopy, in particular, MAC Mode® AFM developed by Keysight Technologies, Inc., entire maps of specific binding sites are obtained by scanning the tethered sensor molecule across the surface to detect binding events (2).

More recently, a Keysight imaging solution known as PicoTREC has been used to simultaneously record two separate images, one that provides the topography of target molecules on a surface and another that displays a map of specifically recognized target molecules on a surface (3, 4). Using MAC Mode, PicoTREC, and optimized AFM cantilevers, a ligand can be kept in close proximity to the surface, allowing efficient recognition and gentle interaction between tip and sample during scanning. PicoTREC resolves molecular recognition during the lateral scan by processing the asymmetric reduction of the oscillation amplitude. In this way, the locations of target molecules are easily determined from their coordinates on the recognition image.

PicoTREC has been demonstrated with an antihistone antibody tethered to the AFM tip recognizing chromatin molecules bound to a mica surface (Figure 2). The recognition image disappears when avidin-biotin interactions are blocked by free biotin (5). This method eliminates requirements for secondary reporter systems. For example, no fluorescence, radioactivity, or enzyme-linked detection schemes are needed when using PicoTREC. Furthermore, the system is extremely sensitive, since single molecule interactions are readily detected. PicoTREC can be used for nanometer-scale epitope mapping of biomolecules, to localize receptor sites during biological processes, or to expedite force spectroscopy studies on areas of interest.

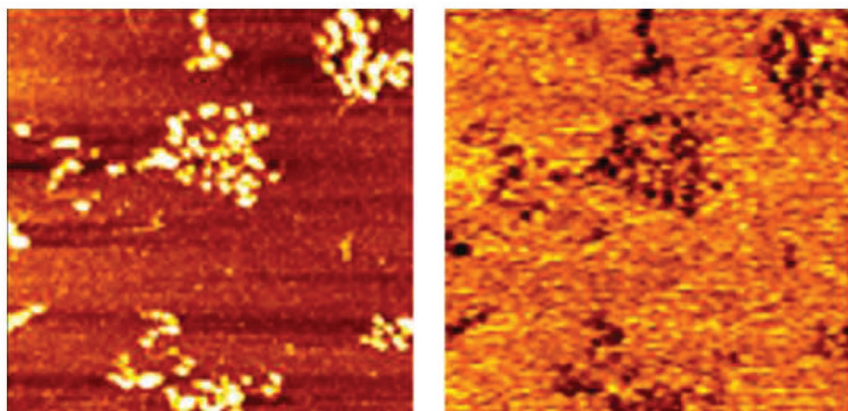


Figure 2. PicoTREC image of MMTV chromatin (DNA and histone protein complex) covalently attached to mica. The Topography image (left) shows individual chromatin molecules. The RECOgnition image (right) shows single molecule binding events between the antihistone antibody, which has been attached to the AFM tip, and histone molecules on the surface. Scan size approximately 500 nm x 500 nm.

## References

1. Hinterdorfer, P., Meth. Cell Biol. **68** (2002) 115-139.
2. Raab, A., Han, W., Badt, D., Smith-Gill, S.J., Lindsay, S.M., Schindler, H., Hinterdorfer, P., Nature Biotech **17** (1999) 902-905.
3. Stroh et al PNAS August 24, 2004 vol. 101 no.34 12503-12507
4. Riener, C.K., Stroh, C.M., Ebner, A., Klampfl, C., Gall, A.A., Romanin, C., Lyubchenko, Y., hinterdorfer, P., Gruber, H.J., Anal. Chim. Acta. **479** (2003) 59-75.

*This application note was created by Cordula M. Stroh, Christian K. Riener, Andreas Ebner, Ferry Kienberger, Hermann J. Gruber, and Peter Hinterdorfer from Johannes Kepler Institute's Institute for Biophysics in collaboration with Keysight Technologies.*

## AFM Instrumentation from Keysight Technologies

Keysight Technologies offers high-precision, modular AFM solutions for research, industry, and education. Exceptional worldwide support is provided by experienced application scientists and technical service personnel. Keysight's leading-edge R&D laboratories are dedicated to the timely introduction and optimization of innovative and easy-to-use AFM technologies.

[www.keysight.com/find/AFM](http://www.keysight.com/find/AFM)

For more information on Keysight Technologies' products, applications or services, please contact your local Keysight office. The complete list is available at: [www.keysight.com/find/contactus](http://www.keysight.com/find/contactus)

### Americas

Canada	(877) 894 4414
Brazil	55 11 3351 7010
Mexico	001 800 254 2440
United States	(800) 829 4444

### Asia Pacific

Australia	1 800 629 485
China	800 810 0189
Hong Kong	800 938 693
India	1 800 112 929
Japan	0120 (421) 345
Korea	080 769 0800
Malaysia	1 800 888 848
Singapore	1 800 375 8100
Taiwan	0800 047 866
Other AP Countries	(65) 6375 8100

### Europe & Middle East

Austria	0800 001122
Belgium	0800 58580
Finland	0800 523252
France	0805 980333
Germany	0800 6270999
Ireland	1800 832700
Israel	1 809 343051
Italy	800 599100
Luxembourg	+32 800 58580
Netherlands	0800 0233200
Russia	8800 5009286
Spain	800 000154
Sweden	0200 882255
Switzerland	0800 805353
	Opt. 1 (DE)
	Opt. 2 (FR)
	Opt. 3 (IT)
United Kingdom	0800 0260637

For other unlisted countries:  
[www.keysight.com/find/contactus](http://www.keysight.com/find/contactus)  
 (BP-09-23-14)

This information is subject to change without notice.  
 © Keysight Technologies, 2007 - 2014  
 Published in USA, July 31, 2014  
 5989-6378EN  
[www.keysight.com](http://www.keysight.com)